

Table Top TDS / Conductivity Meter

ABT202



Product Manual / Warranty Card

Introduction

We thank you for having purchased bench top multi- parameters meter. Before using the instrument, please note that the operation instructions should be read carefully, which will help you to operate and maintain the instrument.


Specification:

Model	ABT202 - Bench Top TDS / Conductivity Meter			
Parameter	Conductivity	TDS	Salt	Temp
Range	0 to 100 mS/cm	0 to 66000 PPM	0 to 50 ppt	0 to 110 Deg C
Accuracy	1%	1%	1%	±0.2
Resolution	1µS / 1ms	1 PPM/PPT	1 PPM	0.1 Deg C
Calibration	0 µS/cm , 84 µS/cm , 1413 µS/cm , 12.88 mS/cm, 80 mS/cm			
ATC	Yes (100 Deg C)			
Battery	9V Battery <i>(Not Supplied with Meter)</i>			
Sensor	Conductivity & Temperature sensors included.			

Accessories:



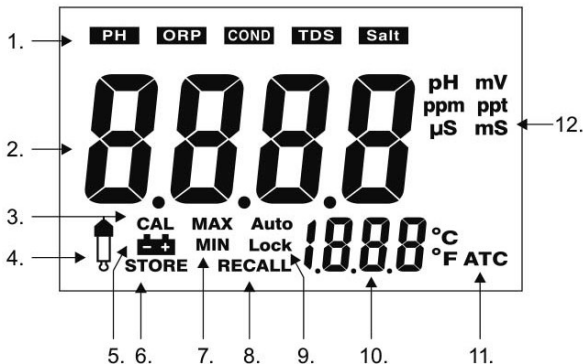


Conductivity Sensor, Temp. Sensor, Standard Solutions,
Power Adapter, Electrode Holder, USB Software,

Display Details



1. Function Mode
2. Measuring Value
3. Calibration Mode
4. Calibration error Indicator
5. Battery Power Low Indicator
6. Reading stored indicator
7. MAX & MIN Value
8. Recall Mode
9. Auto lock or manually lock current reading
10. Temperature Indicator
11. Auto Temperature Compensation
12. Unit

Keyboard Details

	Single Press Lock the current reading, Press for 3 sec. to enter or exit MAX/MIN mode.
	Single Press to Store the current reading. Press 3 sec. to enter Recall mode.
	In Recall mode, browse records. Press both together 3 sec. to enter Advance Setting
	Single Press to Choose different Function mode. Press 3 sec. to switch °C/°F, or switch or switch mg/L or ppm
	Turn on or off power. Press 3 sec. to enter calibration mode

Meter Installation

- 1) Connect AC adaptor with power source.
- 2) Connect the Conductivity Sensor and Temperature Electrode
- 3) Assemble the Electrode Stand and Holder.
- 4) Clean the Electrodes with DM Water and Power On the Meter

Conductivity Calibration:

Switch the Meter to Conductivity Mode. Now Dip the Conductivity electrode and the temperature probe into the 1413 $\mu\text{S}/\text{cm}$ buffer solution and Stir gently until the reading is stable. Press and hold CAL for 3 seconds & the display will appear CAL and flashing 1413. Then the display indicates "SA", then "End" i.e. calibration is done. Meter will return to Measurement Mode.

Note:

1. Calibrated by 12.88 mS/cm standard solution is recommended for measuring high conductivity solution.
2. Calibration error indicator icon will appear, and "Err" instead of "SA", if calibration fails.
3. If the reading is not 0 $\mu\text{S}/\text{cm}$ while the cell is in the air and not dipped into any solution, calibrate it in the air to make reading becomes 0 $\mu\text{S}/\text{cm}$.
4. The other calibration points are 0, 84 $\mu\text{S}/\text{cm}$, 1413 $\mu\text{S}/\text{cm}$, 12.88 mS/cm and 80.0 mS/cm .

Conductivity / TDS / Salt Measurement:

Press **MODE** to choose respective modes.

After calibration, rinse the Conductivity electrode with clean water and wipe it dry. Dip the electrode and the temperature probe into the sample solution which is going to be measured. Stir gently and wait until a stable reading can be obtained.

Note:

- 1.The display will appear “----” when it is over measuring range.
- 2.The unit will auto-range to $\mu\text{S}/\text{cm}$ or mS/cm , or ppm or ppt.
- 3.Don't touch or wipe the surface of the inner black plate of the Sensor
- 4.After measurement, rinse the electrode with clean water. Replace the soaking bottle. The soaking bottle should be always filled with soaking solution

MAX/MIN mode:

Press and hold **MAX/MIN** button until the display appears flashing MAX and MIN icons to enter MAX/MIN. Press **MAX/MIN** lightly to browse MAX and MIN value during this mode. To exit this mode, press and hold **MAX/MIN** button again until the flashing MAX and MIN icons disappear and return to measuring mode.

Store and Recall mode:

In measuring mode, press **STO** to store the current reading. The Store icon and the ordinal of this record will appear on the display.

Press and hold **RCL** for 3 sec. to enter Recall mode. In this mode, use ▲ or ▼ to browse records. Press and hold **STO** to exit this mode and return to measuring mode.

In Recall mode, press ▲▼ together for 3 sec. to clean all the records in the memory.

Advanced Setting:

In TDS mode, press ▲▼ together for 3 sec. to enter advanced setting.

Press **STO** to the ratio between Conductivity and TDS.

Use ▲▼ to adjust ratio (Range 0.5 to 1.0)

Then press **MODE** to confirm and return to measuring mode.

Press **LOCK** to turn on or off “Auto Lock” function. Use ▲▼ to choose.

Then press **MODE** to confirm and return to measuring mode.

In advanced settings, press **CAL** to reset all settings to factory setting.

WARRANTY CARD

Customer Details:

Name: _____

Address: _____

Email Id: _____

Contact No: _____

Dealer Details:

Name: _____

Stamp & Signature

Product Details:

Bill No: _____ Bill Date: _____

Product Serial No: _____

Manufactured By



Rakiro Biotech Systems Pvt Ltd

www.rakiro.net