



BENCH TOP PH / ORP METER

ASM3000



Product Manual / Warranty Card

Introduction

We thank you for having purchased Magnetic Stirrer.

Refer the Product Manual, Datasheet and MSDS before operation.

Specification:

- Model : ASM3000
- Speed Range : 200 – 2000 RPM
- Speed Control : Analog
- Capacity : Up to 3 Liters (H₂O)
- Motor Type : Brushless DC Motor
- Plate Size : 125 x 125 mm
- Power Consumption : 5W – 10W
- Power Supply : 100 – 240V AC, 50/60 Hz
- Power Adapter: Provided 9V1A
- Weight :500 GMS
- Ambient Conditions: 10 – 40°C, ≤80% RH

Accessories:

One Magnetic Stirrer, One Adapter , One Needle

Installation

Place the stirrer on a flat, stable, and level surface.

Connect the power adapter to the stirrer and plug it into the power supply.

Place the stir bar inside the container (beaker/flask).

Position the container on the center of the stirrer.

Note: Ensure the container is placed exactly at the center of the stirrer for smooth and efficient stirring.

Operation

Confirm that installation is completed as per the above instructions.

Turn on the main power supply and switch on the stirrer from the rear side.

Rotate the control knob to start stirring and adjust the speed as required.

After use, turn the knob to the OFF position and switch off the instrument.

Cleaning & Maintenance

Ensure the instrument is switched off and disconnected from the power supply before cleaning.

Clean the outer surface using a soft cloth dampened with water and mild detergent only.

Wipe any spills immediately to prevent damage.

Do not immerse the instrument in water or any liquid.

WARRANTY CARD

Customer Details:

Name: _____

Address: _____

Email Id: _____

Contact No: _____

Dealer Details:

Name: _____

Stamp & Signature

Product Details:

Bill No: _____ Bill Date: _____

Product Serial No: _____

Manufactured By



Rakiro Biotech Systems Pvt Ltd

www.rakiro.net