



RAKIRO BIOTECH SYSTEMS PVT LTD

R-466, TTC Industrial Area, MIDC

Rabale, Navi Mumbai - 400701

Tel No. +91-022-47804040

Email :- sales@rakiro.net

Product Code

BSV1000

Product Name

BACTASLYDE - H2S TEST VIAL (1 Test)

PRODUCT MANUAL

Water contaminated with bacteria is a major cause of water-borne diseases like diarrhoea, dysentery, typhoid, cholera, etc. Detecting these dangerous but unseen predators is quite a tough task and normally it is the domain of the white coated microbiologist who works amidst sterile conditions and a lot of hush hush secrecy. No longer! Just do it yourself with BACTASLYDE. Our unique Water Potability Test Kit detects harmful bacteria with ease and convenience.

Instructions to Use

Step One

Wash hands thoroughly with soap & water before handling the product



Step Two

Remove the Cap gently and pour the water sample up to the mark (20mL)



Step Three

Close the Cap tightly and shake gently to mix properly.

Step Four

Keep the Bottle at Room temperature (20 - 30 Deg C).(No Need of Incubator)

Step Five

Observe the Bottle 24 - 48 Hrs.



**If colour does not change,
the Water is Potable**



**If colour changes to black
the Water is Not Potable**

Note: Conclude the result in 48 hours

Disclaimer: The presentation of H2S Test Kit is based on best of the scientific knowledge published in journals and research papers. It is assumed that the referee and the user is aware of its limitation. Manufacturer does not warrant or assume any type of liability or responsibility for the use & results of this test. Refer Product Literature & MSDS for more details. Single Use Test only. Do not Reuse.